

Distribuido por:



García Martín 19  
28224 Pozuelo, Madrid  
T: 917 091 540 | ventas@anade.com  
anade.com



## ClearKlens Plus

Sterile concentrated neutral detergent for manual use

# VH5

### Description

**ClearKlens Plus** is a concentrated, neutral, high-foaming liquid detergent for the Pharmaceutical industry.

### Key properties

**ClearKlens Plus** is recommended for manual cleaning of process equipment, floors and walls.

**ClearKlens Plus** contains a concentrated blend of high foaming detergents and emulsifiers. Its balanced neutral formulation is particularly effective against product residues found in the Pharmaceutical and Cosmetic industries. It can be used on most types of surfaces including plastics and soft metals such as aluminium.

**ClearKlens Plus** is packaged in 30ml double bagged unit doses, which have been filtered to 0.2µm and sterile filled.

### Benefits

- 0.2µm filtered and gamma irradiated
- Double bagged for transfer into cleanroom
- Suitable for manual use
- Suitable for soft or hard water
- Low level residue detection method available

### Use instructions

Empty the contents of a 30ml bottle of **ClearKlens Plus** into 4 - 10 litres of water

### Technical data

Appearance clear, pale brown liquid  
Relative Density at 20°C 1.03  
pH (1% solution at 20°C) 5.9  
Chemical Oxygen Demand (COD) 400 gO<sub>2</sub>/kg  
Nitrogen Content (N) none  
Phosphorous Content (P) none  
The above data is typical of normal production and should not be taken as a specification.

### Safe handling and storage information

Store in original closed containers or (where applicable) in approved bulk tanks, away from sunlight and extremes of temperature. Full guidance on the handling and disposal of this product is provided in a separate Material Safety Data Sheet.

### Product compatibility

**ClearKlens Plus** is safe for use on materials commonly found in the Pharmaceutical and Cosmetic industries when applied under the recommended conditions. In the event of uncertainty it is advisable to evaluate individual materials before any prolonged use.



## ClearKlens Plus

### Test method

Validated Titration Test method.  
HPLC Test Method.

### Available pack sizes

Article code	Pack size	Pack type
7515165	100x30ml	

### Diversey Europe Operations B.V.

De Boelelaan 32  
1083 HJ Amsterdam  
The Netherlands  
Tel +31 (0)20 258 9700  
[www.diversey.com](http://www.diversey.com)

P92668XFEN-02

© Copyright 2012 Diversey

Distribuido por:



García Martín 19  
28224 Pozuelo, Madrid  
T: 917 091 540 | [ventas@anade.com](mailto:ventas@anade.com)  
[anade.com](http://anade.com)